CAT

Ingenieurbüro CAT
M. Zipperer GmbH

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Sandbath H 4
Sandbad H 4

Table of Content

1	User Instructions.....	4
1.1	Important Instructions for your safety.....	4
2	Danger symbols in the operating manual.....	4
2.1	Danger symbols:	4
2.2	Danger levels	5
2.3	Warning signs on the device.....	5
3	General safety warnings and instructions.....	5
4	General Information	7
5	Scope of Delivery.....	8
6	Intended Use	8
7	Setting up and Starting up the Instrument.....	8
7.1	Unpacking the Instrument.....	8
7.2	Setting up the Instrument	8
7.3	Electrical Connection.....	9
7.4	Switching the instrument on	10
8	Cleaning and Maintenance	11
9	Dismantling, Transport and Storage.....	11
9.1	Dismantling.....	11
9.2	Transport and Storage.....	11
10	Warranty and Liability	12
11	Technical Data.....	12
12	Repairs	12
	Konformitätserklärung / Declaration of conformity.....	20

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1 User Instructions

1.1 Important Instructions for your safety

- Every user must read and understand this manual completely before use. Failure to do so can result in serious injury or death.
- Comply with all safety and accident-prevention regulations applicable to laboratory work.
- Follow general instructions for hazard prevention and general safety instructions, e.g. wear protection clothing, eye protection and gloves.
- This operating manual is part of the product. Thus, it must always be easily accessible.
- This instruction sheet does not purport to address all of the safety problems which might result from the use of this device, chemicals, reagents, apparatus or equipment employed in any specific test or protocols. It is the responsibility of the user to consult their authorized safety advisors and establish appropriate health and safety practices and then determine the application of regulatory limitations prior to use.
- Enclose this operating manual when transferring the device to another place.
- If this manual is lost, please request another one. Please contact your dealer or

Ingenieurbüro CAT
M. Zipperer GmbH
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen
Tel.: ++49-(0)7634-5056-800
Fax: ++49-(0)7634-5056-801
www.cat-ing.de
info@cat-ing.de

2 Danger symbols in the operating manual

The safety instructions in this manual appear with the following danger symbols and danger levels:

2.1 Danger symbols:

 Hazard point	 Electrical shock
 Risk of fire	 Explosion
 Bio hazard	 Hot surface
 Chemical hazard	 Material damage

2.2 Danger levels

 DANGER	Will lead to severe injuries or death
 WARNING	May lead to severe injuries or death
 CAUTION	May lead to light to moderate injuries
 NOTICE	May lead to material damage

2.3 Warning signs on the device

WARNING

This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

WARNING

This symbol indicates that there are **hot surfaces** on the device. It refers to the hotplate of the device which can reach up to 500°C during operation causing a hazard resulting in serious burns.

- Avoid contact with the hotplate, vessel and heating medium.
- Avoid storage of flammable substances near the instrument.

3 General safety warnings and instructions

	<p> DANGER Risk of explosion.</p> <ul style="list-style-type: none"> • Do not operate the device in the vicinity of highly flammable or explosive substances. The instrument is not explosion-proof. • Do not use this device for processing any substances which could generate an explosive atmosphere. • Do not use this device to process any explosive or highly reactive substances. • Do not use flammable or explosive substances near the instrument.
	<p> DANGER Electric shock as a result of penetration of liquid.</p> <ul style="list-style-type: none"> • Do not allow any liquids to penetrate the inside of the housing • Switch off the device and disconnect the power plug before starting cleaning or disinfection work. The On/Off Switch on the device does not disconnect the device from the power source. • Use only closed tubes. • Only plug the device back in if it is completely dry, both inside and outside.

	<p>▲WARNING Risk from incorrect supply voltage</p> <ul style="list-style-type: none"> ● Only connect the device to an AC power source with an protective earth (PE). ● Only connect the device to voltage sources which correspondent to the electrical requirements on the type label.
	<p>▲WARNING Electric shock due to damage to device or mains cable</p> <ul style="list-style-type: none"> ● Only connect the device to the mains supply if the device and the mains cable are undamaged ● Only use devices that have been properly installed or repaired. ● In case of danger, disconnect the device from the mains supply by pulling the power plug from the mains socket or by using the isolating device intended for this purpose (e.g. emergency stop switch)
	<p>▲WARNING Lethal voltage inside the device</p> <ul style="list-style-type: none"> ● Do not open the device. ● Ensure that the housing is always closed and undamaged so that no parts inside the housing can be contacted by accident. ● The On/Off Switch on the device does not disconnect the device from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely. ● Do not allow any liquids to penetrate the inside of the housing. ● Repairs are only to be carried out by trained service technicians.
	<p>▲WARNING Damage to health due to corrosive or aggressive chemicals</p> <ul style="list-style-type: none"> ● Observe all markings on the reagent bottles. ● Always check the instrument for leaks and air bubbles. Special attention should be directed to determine that all push-ons, threaded connections and suction tubes are firmly in place before beginning operation. Leaking solutions may endanger persons and materials ● Use proper connecting vessels, protective clothing and gloves. ● Avoid splashes ● When dispensing, maintain a physical distance between the instrument and the body. ● Dangerous and fuming chemicals must be dispensed in a fume hood. ● Only employ the instrument for the purpose intended by the manufacturer, and particularly within the resistance limits of the instrument. If in doubt, contact your supplier, or the manufacturer's factory representative at the phone number shown at the front page of this operating instruction. ● Always use the instrument in such a manner that neither the operator, nor any other person is endangered.
	<p>▲WARNING Damages to health due to infectious liquids and pathogenic germs.</p> <ul style="list-style-type: none"> ● When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes. ● Wear personal protective equipment ● For comprehensive regulations about handling germs or biological material of the risk group II or higher, please refer to the "Laboratory Biosafety Manual" in its respectively current valid version from the World Health Organisation

	<p>▲WARNING Damage to health due to contaminated device and accessories</p> <p>In the following cases, sample material can be released:</p> <ul style="list-style-type: none"> - improperly sealed tubes - unstable tubes - high vapour pressure of the content so that the seal of the tubes can spring open - damaged sealings - smashed glass tubes <ul style="list-style-type: none"> ● Observe the nationally prescribed safety environment when working with hazardous, toxic and pathogenic samples. Pay particular attention to personal protective equipment (gloves, clothing, goggles, etc.), extraction, and the safety class of the lab. ● Decontaminate the device and the accessories before storage and shipping.
	<p>▲WARNING Risk of fire</p> <ul style="list-style-type: none"> ● Do not use this device to process any highly flammable liquids
	<p>▲CAUTION Poor safety due to inadequate fixing of the unit</p> <ul style="list-style-type: none"> ● Ensure that the unit is firmly attached to a solid stand.
	<p>▲CAUTION Poor safety due to incorrect accessories and spare parts.</p> <p>The use of accessories and spare parts other than recommended by Ingenieurbüro CAT, M. Zipperer GmbH may impair the safety, function and precision of the device. Ingenieurbüro CAT, M. Zipperer GmbH cannot be held liable or accept any liability for damage resulting from the use of incorrect or non-recommended accessories and spare parts, or from the improper use of such equipment.</p> <ul style="list-style-type: none"> ● Only use accessories and spare parts recommended by Ingenieurbüro CAT, M. Zipperer GmbH

4 General Information

Our Hotplates are designed according to protection class 1. They were manufactured and tested according to DIN EN 61010. According to these regulations, the devices are designed to meet the requirements for safe and correct operations. To ensure the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

▲WARNING It is the user's task to find out whether the device is suitable for the application. If in doubt clarify this with your dealer or directly with the manufacturer.

▲WARNING Please comply with all safety and accident-prevention regulations applicable to laboratory work.

5 Scope of Delivery

Inspect the content of the package for damage or missing parts:

1 Heater H 4 (Part no.: 60224-0000)
1 Instruction Manual

6 Intended Use

This hotplate is mainly to be used to heat hydrous solutions and oil not containing inflammable parts in glass vessels. It is designed for use in chemical and biological laboratories of industrial enterprises, universities and pharmacies.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

The onus is on the user to find out whether the device is suitable for his application. If in doubt clarify this with your dealer or with the manufacturer directly.

7 Setting up and Starting up the Instrument

7.1 Unpacking the Instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

If the instrument is not damaged and all parts are complete you may start to operate the device after reading the instruction manual.

7.2 Setting up the Instrument

Please place the device on a fire-proof horizontal and even surface. A minimum distance of 50cm to inflammable materials should always be observed.

▲ DANGER The device may not be operated in explosion-prone areas.

▲ WARNING The power cable may not come in contact with the hotplate.

▲ WARNING The device is not to be used without supervision.

▲ WARNING When connecting the instrument to an AC power outlet ensure that your local supply voltage matches the indication on the instrument.

7.3 Electrical Connection

▲WARNING 

- The units are equipped with a grounded power cord of protection class 1, IEC 884-1 (DIN 49441 R2) 16 A, 250 V. For North America, with a U.S. standard plug (NEMA 5 ,Type B) 10 A, 125 V.
- When operating the instruments in countries with different AC plug systems use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.
- The instrument is earthed as supplied. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug!
- When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.
-

Before connection the instrument to the mains ensure that the rocker switch (1) is set to the „OFF“- position.

▲WARNING

When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.

7.4 Switching the instrument on

- Start heating with rocker switch (1)
- Set desired heating performance with the temperature control button (2).
- The temperature is continuously adjustable from 0 – 330°C
- "Ready for use" is indicated by the left signal lamp (3) as soon as the heating plate is supplied by power.

▲WARNING **Hot!** The device has hot areas especially the stainless steel bath itself which can reach a temperature of 330°C. Hazard of serious burns. Avoid contact with bath container, sand and liquid. Also avoid flammable substances to be stored near the instrument.

▲WARNING The sandbath will stay hot for about 15 minutes after being switched off. Hazard of serious burns. Avoid contact with hotplate, vessel and liquid. Also avoid flammable substances to be stored near the instrument.

Notes on setting the hotplate temperature:

The contents of the vessel will always have a lower temperature than the set hotplate temperature. This is a result of the differences in heat conductivity of various vessels, volume of medium to be heated and heat loss due to radiation.

The hotplate temperature therefore should be set 1.5 to 4 times that of the desired temperature of the medium. The value should be increased according to the specific heat capacity, volume and desired temperature of the medium.

8 Cleaning and Maintenance

The casing and the stainless steel container of the sandbath are mostly acid resistant. The heating surface stays constantly smooth and is easy to clean. Slightly soiled surfaces can be cleaned (when hand warm or cold) with water and dish washing liquid. Do not use steel wool or a similar scratching sponge to clean the units.

Crust, limestone and water stains may be cleaned with the usual household detergents. Please remove remaining detergent thoroughly with a damp cloth or sponge, since some of the detergents react caustic with higher temperatures. If soils are already encrusted use a razorblade scratcher. Plastics, kitchen foil as well as stuff containing sugar must be removed immediately.

Mechanical supports are: Razorblade scratchers, non-scratching sponges and non-scratching steel wool.

The device is maintenance-free. Should repairs become necessary, only trained and authorized service personnel may open and repair the instrument.

Unplug the AC power cord before opening the housing.

9 Dismantling, Transport and Storage

9.1 Dismantling

1. Switch off the unit by using switching the rocker switch (1) to the „OFF“-Position.
2. Disconnect the unit from the mains.
3. Now you may remove the instrument from the working area.

9.2 Transport and Storage

Prior to transport:

Switch the instrument off and unplug the power supply.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, please sort according to materials: metal, glass, plastic, etc. Also be sure to dispose of the packing material in an environmental-friendly manner.

10 Warranty and Liability

The manufacturer agrees to correct for the original user of this product, either by repair, or at the manufacturer's election, by replacement, any defects in material or workmanship which develop within 24 months after delivery of this product to the original user. In the event of replacement, the replacement unit will be warranted for the remainder of the original twenty-four (24) month period of ninety (90) days, whichever is longer.

If this product should require service, contact your distributor or manufacturer for necessary instructions.

This warranty shall not apply if the defect or malfunction was caused by accident, neglect, unreasonable use, improper service, or other causes not arising out of defects in material or workmanship.

There are no warranties, expressed or implied, including, but not limited to, those of merchantability or fitness for a particular purpose, which extended beyond the description and period set forth herein.

The manufacturer's sole obligation under this warranty is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind resulting from use or possession of the product.

11 Technical Data

Model	Part. No.	Vessel Size (mm)	Material	Heating Power in Watt	Temperature Range in °C (Hotplate)	Dimension (WxDxH) in mm	Weight (Kg)
H 4	60224-0000	140x140x55	Stainless steel	600	40-330	155x157x165	1,8

Power Supply : see rating plate on device

Condition of Surrounding:

Permissible ambient temperature : +5 °C to + 40 °C
Permissible ambient temperature : max 80% relative humidity

12 Repairs

WARNING

When returning instruments for repair that have come into contact with hazardous substances, please

Provide precise information on the relevant medium

Take protective measures to ensure the safety of our receiving and maintenance personnel

Mark the package as appropriate for hazardous materials.

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