CAT

Ingenieurbüro CAT
M. Zipperer GmbH

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Magnetic Stirrer
Magnetrührer
ECM 2, ECM 5, M 5, M 15

Content

1	User Instructions	4
1.1	Important Instructions for your safety	4
1.2	Danger symbols in this operating manual	4
1.2.1	Danger symbols:.....	4
1.2.2	Danger levels.....	5
1.2.3	Warning signs on the device	5
2	General safety warnings and instructions	6
3	General Information.....	8
4	Intended Use	8
5	Setting up and Starting up the Instrument.....	9
5.1	Unpacking the Instrument	9
5.2	Setting up the Instrument.....	9
6	Electrical Connection.....	10
6.1	Switching the instrument on.....	10
7	Cleaning and Maintenance.....	11
8	Dismantling, Transport and Storage.....	11
8.1	Dismantling	11
8.2	Transport and Storage	11
9	Warranty and Liability	12
10	Technical Data	13
11	Repairs.....	13
12	Repair Return Form	14
25	Declaration of Conformity / Konformitätserklärung.....	24

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1 User Instructions

1.1 Important Instructions for your safety

- Every user must read and understand this manual completely before use. Failure to do so can result in serious injury or death.
- Comply with all safety and accident-prevention regulations applicable to laboratory work.
- Follow general instructions for hazard prevention and general safety instructions, e.g. wear protection clothing, eye protection and gloves.
- This operating manual is part of the product. Thus, it must always be easily accessible.
- This instruction sheet does not purport to address all of the safety problems which might result from the use of this device, chemicals, reagents, apparatus or equipment employed in any specific test or protocols. It is the responsibility of the user to consult their authorized safety advisors and establish appropriate health and safety practices and then determine the application of regulatory limitations prior to use.
- Enclose this operating manual when transferring the device to another place.
- If this manual is lost, please request another one. Please contact your dealer or

Ingenieurbüro CAT
M. Zipperer GmbH
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen
Tel.: ++49-(0)7634-5056-800
Fax: ++49-(0)7634-5056-801
www.cat-ing.de
info@cat-ing.de

1.2 Danger symbols in this operating manual

The safety instructions in this manual appear with the following danger symbols and danger levels:

1.2.1 Danger symbols:

	Hazard point		Electrical shock
	Risk of fire		Explosion
	Crushing		Hot surface

	Bio hazard		Magnetic field
	Chemical hazard		Material damage

1.2.2 Danger levels

	Will lead to severe injuries or death
	May lead to severe injuries or death
	May lead to light to moderate injuries
	May lead to material damage

1.2.3 Warning signs on the device

WARNING

This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

WARNING

Beware of the effect of the **magnetic field** on cardiac pacemakers and data storage media.

WARNING

This symbol indicates that there are **hot surfaces** on the device. It refers to the hotplate of the device which can reach up to 500°C during operation causing a hazard resulting in serious burns. • Avoid contact with the hotplate, vessel and heating medium.

- Avoid storage of flammable substances near the instrument.

2 General safety warnings and instructions

	<p>⚠ DANGER Risk of explosion.</p> <ul style="list-style-type: none"> • Do not operate the device in the vicinity of highly flammable or explosive substances. The instrument is not explosion-proof. • Do not use this device for processing any substances which could generate an explosive atmosphere. • Do not use this device to process any explosive or highly reactive substances. • Do not use flammable or explosive substances near the instrument.
	<p>⚠ DANGER Electric shock as a result of penetration of liquid.</p> <ul style="list-style-type: none"> • Do not allow any liquids to penetrate the inside of the housing • Switch off the device and disconnect the power plug before starting cleaning or disinfection work. The On/Off Switch on the device does not disconnect the device from the power source. • Use only closed tubes. • Only plug the device back in if it is completely dry, both inside and outside.
	<p>⚠ WARNING Risk from incorrect supply voltage</p> <ul style="list-style-type: none"> • Only connect the device to a AC power source with an protective earth (PE). • Only connect the device to voltage sources which correspondent to the electrical requirements on the type label.
	<p>⚠ WARNING Electric shock due to damage to device or mains cable</p> <ul style="list-style-type: none"> • Only connect the device to the mains supply if the device and the mains cable are undamaged • Only use devices that have been properly installed or repaired. • In case of danger, disconnect the device from the mains supply by pulling the power plug from the mains socket or by using the isolating device intended for this purpose (e.g. emergency stop switch)
	<p>⚠ WARNING Lethal voltage inside the device</p> <ul style="list-style-type: none"> • Do not open the device. • Ensure that the housing is always closed and undamaged so that no parts inside the housing can be contacted by accident. • The On/Off Switch on the device does not disconnect the device from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely. • Do not allow any liquids to penetrate the inside of the housing. • Repairs are only to be carried out by trained service technicians.
	<p>⚠ WARNING Damage to health due to corrosive or aggressive chemicals</p> <ul style="list-style-type: none"> • Observe all markings on the reagent bottles. • Always check the instrument for leaks and air bubbles. Special attention should be directed to determine that all push-ons, threaded connections and suction tubes are firmly in place before beginning operation. Leaking solutions may endanger persons and materials

	<ul style="list-style-type: none"> ● Use proper connecting vessels, protective clothing and gloves. ● Avoid splashes ● When dispensing, maintain a physical distance between the instrument and the body. ● Dangerous and fuming chemicals must be dispensed in a fume hood. ● Only employ the instrument for the purpose intended by the manufacturer, and particularly within the resistance limits of the instrument. If in doubt, contact your supplier, or the manufacturer's factory representative at the phone number shown at the front page of this operating instruction. ● Always use the instrument in such a manner that neither the operator, nor any other person is endangered.
	<p>▲WARNING Damages to health due to infectious liquids and pathogenic germs.</p> <ul style="list-style-type: none"> ● When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes. ● Wear personal protective equipment ● For comprehensive regulations about handling germs or biological material of the risk group II or higher, please refer to the "Laboratory Biosafety Manual" in its respectively current valid version from the World Health Organisation
	<p>▲WARNING Damage to health due to contaminated device and accessories</p> <p>In the following cases, sample material can be released:</p> <ul style="list-style-type: none"> - improperly sealed tubes - unstable tubes - high vapour pressure of the content so that the seal of the tubes can spring open - damaged sealings - smashed glass tubes <ul style="list-style-type: none"> ● Only mix in closed tubes ● Observe the nationally prescribed safety environment when working with hazardous, toxic and pathogenic samples. Pay particular attention to personal protective equipment (gloves, clothing, goggles, etc.), extraction, and the safety class of the lab. ● Decontaminate the device and the accessories before storage and shipping.
	<p>▲WARNING Risk of fire</p> <ul style="list-style-type: none"> ● Do not use this device to process any highly flammable liquids
	<p>▲CAUTION Poor safety due to inadequate fixing of the unit</p> <ul style="list-style-type: none"> ● Ensure that the unit is firmly attached to a solid stand.
	<p>▲CAUTION Poor safety due to incorrect accessories and spare parts.</p> <p>The use of accessories and spare parts other than recommended by Ingenieurbüro CAT, M. Zipperer GmbH may impair the safety, function and precision of the device. Ingenieurbüro CAT, M. Zipperer GmbH cannot be held liable or accept any liability for damage resulting from the use of incorrect or non-recommended accessories and spare parts, or from the improper use of such equipment.</p> <ul style="list-style-type: none"> ● Only use accessories and spare parts recommended by Ingenieurbüro CAT, M. Zipperer GmbH

▲CAUTION Crush hazard due to moving parts

- Do not replace any consumables as long as the device is running.
- Do not open the coverage as long as the device is running

3 General Information

Our Magnetic Hotplate Stirrers are designed according to protection class 1. They were manufactured and tested according to DIN EN 61010. According to these regulations, the devices are designed to meet the requirements for safe and correct operations. To ensure the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

4 Intended Use

The magnetic stirrer is mainly to be used to stir, oil and other hydrous solutions not containing inflammable parts in glass vessels. It is designed for use in chemical and biological laboratories of industrial enterprises, universities and pharmacies.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

▲WARNING It is the user's task to find out whether the device is suitable for the application. If in doubt clarify this with your dealer or directly with the manufacturer.

▲WARNING Please comply with all safety and accident-prevention regulations applicable to laboratory work.

5 Setting up and Starting up the Instrument

5.1 Unpacking the Instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

If the instrument is not damaged and all parts are complete you may start to operate the device after reading the instruction manual.

5.2 Setting up the Instrument

Please place the device on a fire-proof horizontal and even surface. A minimum distance of 50cm to inflammable materials should always be observed.

▲ DANGER

The device may not be operated in explosion-prone areas.

▲ WARNING

The power cable may not come in contact with the hotplate.

▲ WARNING

The device is not to be used without supervision.

▲ WARNING

When connecting the instrument to an AC power outlet ensure that your local supply voltage matches the indication on the instrument.

▲ WARNING

Beware of the effect of the magnetic field on cardiac pacemakers and data storage media.

6 Electrical Connection

⚠ WARNING ⚡

- A EURO AC plug (DIN 49441 CEE 7/VII10/ 16 A 250 V) is standard on all instruments. For North America instruments feature standard US plugs (NEMA Pub.No.WDI1961 ASA C 73.1. 1961 page 8 15A 125V).
- When operating the instruments in countries with different AC plug systems use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.
- The instrument is earthed as supplied. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug!
- When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.

6.1 Switching the instrument on

	<p>⚠ WARNING Before connecting the instrument to the mains ensure that the rocker switch (1) is set to the „OFF“- position and the speed control button is set to “0”</p>
---	--

1. Switch on the unit with rocker switch (1). It is now illuminated.
2. Set desired speed at control knob (2). Always start at a low speed to avoid splashes.
The speed can be set steplessly for

ECM 2:10 - 200 rpm
 ECM 5:50 - 1200 rpm
 M 5: 60 - 1600 rpm
 M 15: 60 - 1600 rpm

speed control knob (2)

rocker switch (1)

Note: The vessel (e.g. a glass beaker) must be permeable to magnetic field lines.

Note: Best fitting length of stirring bar for:

M 5 : 20 - 60 mm
 M 15 : 20 - 60 mm
 ECM 2 : 20 - 30 mm
 ECM 5 : 20 - 30 mm

	<p>⚠ WARNING When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.</p>
---	--

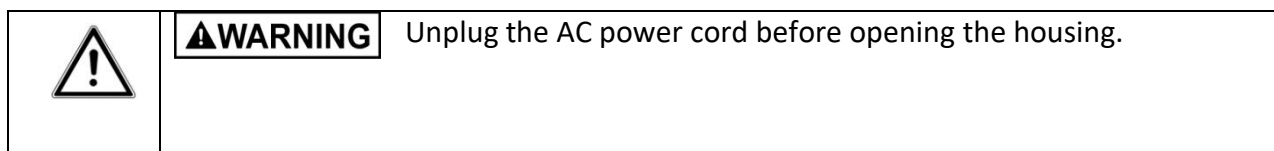
7 Cleaning and Maintenance

The stainless steel casing of the magnetic stirrer ECM 2 (ECM 5, M 5 M 15) is mostly acid resistant. The surface stays constantly smooth and is easy to clean. Slightly soiled surfaces can be cleaned with water and dish washing liquid. Do not use steel wool or a similar scratching sponge to clean the units.

Crust, limestone and water stains may be cleaned with the usual household detergents. Please remove remaining detergent thoroughly with a damp cloth or sponge. If soils are already encrusted use a razorblade scratcher.

Mechanical supports are: Razorblade scratchers, non-scratching sponges and non-scratching steel wool.

The device is maintenance-free. Should repairs become necessary, only trained and authorized service personnel may open and repair the instrument.

8 Dismantling, Transport and Storage

8.1 Dismantling

1. Switch off the unit by switching the rocker switch (1) to the „OFF“-Position.
2. Set speed to “0” position by turning the speed control knob to “0”.
3. Disconnect the unit from the mains.
4. Now you may remove the instrument from the working area.

8.2 Transport and Storage

Prior to transport:

Switch the instrument off and unplug the power supply.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, please sort according to materials: metal, glass, plastic, etc. Also be sure to dispose of the packing material in an environmental-friendly manner.

WEEE No. DE 54398924

9 Warranty and Liability

The manufacturer agrees to correct for the original user of this product, either by repair, or at the manufacturer's election, by replacement, any defects in material or workmanship which develop within 24 months after delivery of this product to the original user. In the event of replacement, the replacement unit will be warranted for the remainder of the original twenty-four (24) month period of ninety (90) days, whichever is longer.

If this product should require service, contact your distributor or manufacturer for necessary instructions.

This warranty shall not apply if the defect or malfunction was caused by accident, neglect, unreasonable use, improper service, or other causes not arising out of defects in material or workmanship.

There are no warranties, expressed or implied, including, but not limited to, those of merchantability or fitness for a particular purpose, which extended beyond the description and period set forth herein.

The manufacturer's sole obligation under this warranty is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind resulting from use or possession of the product.

10 Technical Data

WARNING

The onus is on the user to find out whether the device is suitable for his application. If in doubt clarify this with your dealer or with the manufacturer directly.

Model	Part. No.	Power supply	Speed (rpm)	Motor (W)	Stirring Capacity related to H ₂ O	Working Area (mm)	Dimension (WxDxH) in mm	Weight (Kg)
ECM 2	60252-0000 (60252-0001)	100 – 250 V AC 100 – 250 V AC	10 - 200	6	1 l	100 x 100	100 x 115 x 58	0.7
ECM 5	60255-0000 (60252-0001)	100 – 250 V AC 100 – 250 V AC	50 -1200	6	1 l	100 x 100	100 x 115 x 58	0.7
M 5	60265-0000 60265-0001	230 V AC 115 V AC	60 -1600	25	10 l	150 x 150	157 x 157 x 80	1.8
M 15	60261-0000	230 V AC	60 - 1600	25	10 l	145 x 160	145 x 184 x 85	1,7

Power Supply: see rating plate on device

Condition of Surrounding:

Permissible ambient temperature: +5 °C to + 40 °C

Permissible ambient temperature: max 80% relative humidity

11 Repairs

WARNING

When returning instruments for repair that have come into contact with hazardous substances, please
 Provide precise information on the relevant medium
 Take protective measures to ensure the safety of our receiving and maintenance personnel
 Mark the package as appropriate for hazardous materials.

12 Repair Return Form

Important:

- Calibration Service is automatically performed for instruments that require adjustments
- We contact you by sending a repair estimate.
- Repairs will only be carried out after your release.

CONTACT/USER INFORMATION

Contact:	Phone No.:
Fax No.:	Email:
Billing:	Shipping:
Company	Company
Address	Address

INSTRUMENT INFORMATION

Model	Serial No.
Please describe all problems/malfunctions	

Operating Conditions (please fill in if applicable)

Ambient Temp.	Humidity	Speed
Load	Volume	Viscosity
Temperature in °C	Sample Temperature	Operating Time
Sample Description*		

*NOTE: If the instrument was exposed to hazardous material, it must be decontaminated BEFORE returning it to Ingenieurbüro CAT, M. Zipperer GmbH and an MSDS for hazardous material must be included with the instrument.

RETURN SHIPPING

UPS <input type="checkbox"/>	Air Parcel Post <input type="checkbox"/>	Collect** <input type="checkbox"/>	Other** <input type="checkbox"/>
------------------------------	--	------------------------------------	----------------------------------

**Your account number is required for UPS collect respectively the address and contact of your preferred forwarder if you choose any other transport means.

PACKAGING INSTRUCTIONS TO RETURN A INSTRUMENT FOR REPAIR

- Remove all accessories (e.g. homogeniser tools, stirring paddles, stands, clamps, cables) from the instrument
- Clean excess testing material off the instrument/accessory
- Include MSDS sheets for all hazardous materials used with this instrument
- Pack the instrument in its original box. If the box is not available, take care to wrap the instrument and accessories with enough material to support them.
- DO NOT send accessories unless there is a problem with them. When sending back, wrap each item to avoid contact with the instrument.
- Pack the instrument and related items in a strong box for shipping. Mark the outside of the box with handling instructions.

Example: "Handle with care" or "Fragile- Delicate Instrument" and send to

Ingenieurbüro CAT, M. Zipperer GmbH
 Service Department
 Wettelbrunner Str. 6
 D-79282 Ballrechten-Dottingen

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We declare under our sole responsibility that this product (see table) corresponds to the regulations and conforms with the standards or standardized documents:

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The agreement with further valid guidelines/regulations following for the product is explained:

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Applied one harmonized standards in particular:

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In the case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

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**Ingenieurbüro CAT, M. Zipperer GmbH, Service Department,
Ingenieurbüro CAT M. Zipperer GmbH